Safety Instructions

byko-visc CP Family

Type 300 003 76

Table of Contents

English	4
Deutsch.....	6
Français.....	8
Italiano	10
Español	12
Portuguese	14
Nederlands	16
Dansk	18
Svenska	20
Norsk	22
Suomi	24
Polski	26
Česky	28
Slovenská	30
Slovenščina	32
Magyarul.....	34
România	36
Български	38
ελληνικά	40
Türkçe	42
Русский.....	44
Eesti	46
Latviešu	48
Lietuviškas	50
中文	52
日本語.....	54

Safety Instructions

Before operation read the safety instructions in the instrument manual. No product liability or warranty claims can be made if the device is not operated in accordance with the operating instructions.

If you pass this instrument to somebody else, make sure to include these instructions.

Any questions? Please contact our BYK-Gardner Customer Service:

<https://www.byk-instruments.com>

DANGER: Injuries possible

Defects and extraordinary loads: If safe operation can no longer be presumed, shut down the device and secure it against unintended operation.

The device must be presumed unsafe to operate:

- If visible damage is evident,
- if the instrument is no longer working,
- if it has been stored for long periods under adverse conditions or
- after harsh treatment during shipping.

Do not perform any repairs on the unit. The unit must be opened by BYK-Gardner authorized repair locations only.

In case of operation using the external power supply unit, it is to be ensured the rated voltage for the power supply unit (see rating plate on the power supply unit) matches the mains voltage on your mains socket.

How to disconnect the instrument from any power source:

- By unplugging the mains plug on the mains cable for the external power supply unit or
- by unplugging the DC input from the external power supply unit to the instrument.

Use only the power supply included with delivery.

WARNING: Severe material damage

Only devices meeting the requirements for low voltage safety may be connected to the USB interface.

WARNING! This manual cannot address all of the safety considerations associated with its use. It is the responsibility of the user to consult this manual and establish appropriate safety practices for use with this equipment and the individual material being tested.

WARNING! The **byko-visc** viscometers are designed and intended for the use described in the manual. Using the viscometer for other purposes for which it was not designed may reduce or eliminate the protection offered by the features of the device. Serious injury may result.

WARNING! This instrument is not intended for use in a potentially hazardous environment or with materials that may release toxic or flammable gases at the temperatures encountered during testing.

WARNING! Never defeat the ground conductor or operate the equipment in the absence of a suitably installed ground conductor. Contact the appropriate electrical inspection authority or an electrician if you are uncertain that suitable grounding is available.

WARNING! The sample plate at the bottom of the device heats up during the test procedure and cools down again (*premium models*). Never touch the sample plate or parts that come into contact with the plate during or immediately after the test without appropriate protective equipment. The full measurement range of the instrument is from 5°C to 235°C; the exact range depends on the model used. The temperature of the sample plate is always shown on the display. Before starting any cleaning, wait until the instrument has cooled down to room temperature.

NOTICE

Please note the following points:

- Familiarize yourself with the layout and operation of the controls.
- Ensure that no clothing, jewelry or hair can be caught up in moving parts during use.

Additional information on disposal

This symbol means: Do not dispose of this product together with your household trash. Please refer to the information of your local community or contact our dealers regarding the proper handling of end-of-life electric and electronic equipment. Recycling of this product will help to conserve natural resources and prevent potential negative consequences for the environment and human health caused by inappropriate waste handling.

Technical data

Temperature range	10° C to 40° C (50° F to 104° F) for operation 0° C to 60° C (32° F to 140° F) for storage
Relative humidity	Up to 85% non-condensing at 35° C (95° F)
Operation altitude	Up to 2000 m (6561 ft)
Dimensions (LxWxH)	300 x 170 x 430 mm (11.5 x 6.5 x 17 in)
Weight	6.5 kg (14.3 lbs)
Interface	2x USB Type A 1x USB Type C 1x Ethernet
Power supply instrument	byko-visc CP L models: 24 V  , max. 3.75 A byko-visc CP H models: 24 V  , max. 5 A
External power supply	byko-visc CP L models: Input: 100 - 240 V  ; 50 - 60 Hz ; max. 1.3 A Output: 24 V  ; max. 3.75 A ; max. 90 W byko-visc CP H models: Input: 100 - 240 V  ; 50 - 60 Hz ; max. 1.4 A Output: 24 V  ; max. 5 A ; max. 120 W

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BYK-Gardner GmbH