byko-visc CP

Short Instructions

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Instructions rapides

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Instrucciones resumidas

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Table of Contents

English	4
Deutsch.....	14
Français.....	24
Italiano	34
Español	44
Portuguese	54
中文	64
日本語	74

Table of Contents

1	System Description	5
2	Instrument Setup	6
3	Main Menu	7
4	Measurement Settings	8
5	System Configuration	9
6	Viscosity Measurement	10
7	Browse Measurements	12
8	Technical Data	13

1 System Description

The entire system consists of instrument and power supply. The cone is to be ordered separately and will be available in the packaging.

CAUTION!


Read the safety instructions and the full instruction manual - available via download - before using this instrument. The viscometer has moving parts which should never be touched during operation. Cone and plate can become very hot. Allow to cool down before touching them.

Viscometer Front

- 1 - Housing
- 2 - Touchscreen
- 3 - Moving head
- 4 - Cone attachment
- 5 - Cone (see next page)
- 6 - Plate

Viscometer Back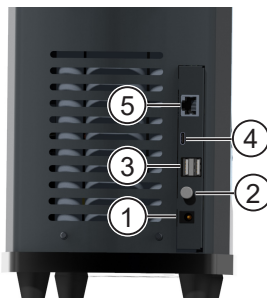

- 1 - External power supply
- 2 - Power button: On and off
- 3 - USB type A:
 - USB flash drive
 - USB printer (or vice versa)
- 4 - USB type C (*)
- 5 - LAN interface (*)

(*) For PC communication (not used yet)


NOTICE

Retain the packaging in case the instrument needs to be shipped. Always drive the viscometer head down before shipping the instrument.

2 Instrument Setup

1. Place the viscometer in a suitable location.
2. Plug the power supply into a mains outlet and plug the output cable into the back of the viscometer head.
3. Turn instrument on with the power button.
4. The viscometer head moves to the upper position.
5. The main menu is shown in the touchscreen display.
6. Hold cone attachment with one hand and insert cone with the other hand.
7. Use the two screws to line up the groove; turn cone clockwise to attach it.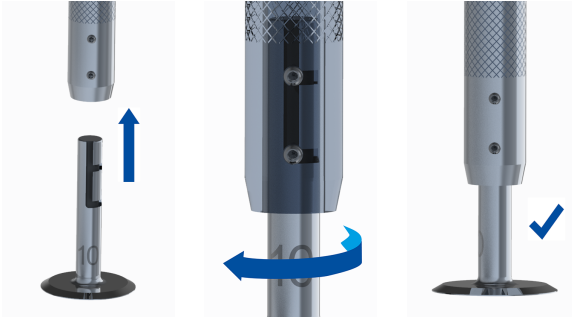

The cone is correctly fastened when one screw fits into each groove.

NOTICE

Check number on the cone and use the **Measurement Settings** to select the cone attached. Use the **Configuration** menu to move the viscometer head up and down, if required.

3 Main Menu

In the main menu the following symbols are displayed.

1 Viscosity Measurement

Perform measurements using existing measurement settings and save results in instrument memory.

2 Browse Measurements

View and manage measurement results saved in instrument memory.

3 Measurement Settings

Change measurement settings and / or instrument configuration.


CAUTION!

Always apply a small amount of sample to the plate before starting a viscosity measurement - i.e. before the cone begins to rotate on the plate.

4 Measurement Settings

Tap the settings icon in the main menu to manage settings changed **often**.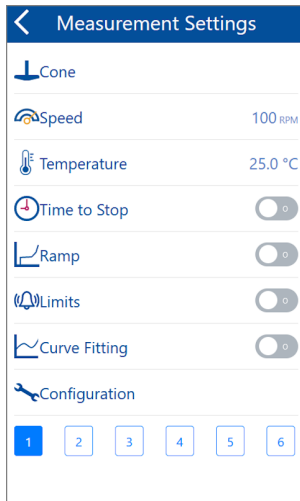

Cone	Select attached cone. Only cones applicable to your model will be shown.
Speed	Input a value from 0.1 - 250 rpm. Cone and speed combination will show the max. viscosity that can be measured with this combination.
Temperature	Set measurement temperature at which the unit will start when reached. Maximum value depends on model.
Time to Stop	Set measurement time in hours / minutes / seconds.
Ramp	Define measurement ramp: Enter start speed, end speed, time to run.
Limits	Select alarm criteria: Set upper limit, set lower limit, select alarm type (buzzer, display flash, both, or none).
Curve Fitting	Choose between: Newton, Bingham, Casson, Ostwald, Hershel-Buckley to overlay onto the measurement graph.
Configuration	Check and change system settings, see System Configuration .
Presets (1 ... 6)	Settings are automatically stored in selected slot. Tap a slot to recall stored settings.

Use back navigation icon (<) to go back to previous page. Current settings will be saved.


NOTICE

Settings for **Speed** and **Cone** will be used as defaults and can also be changed directly during measurements.

5 System Configuration

Tap on the configuration icon in the **Measurement Settings** menu to manage settings changed **seldom**.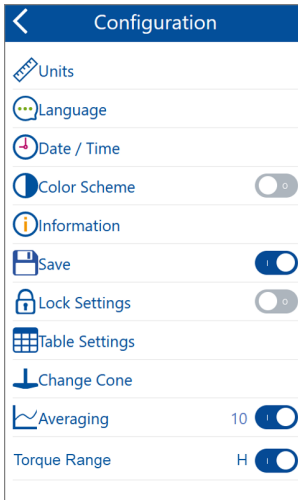

Units	Select units for display of viscosity and temperature.
Language	Select preferred display language.
Date / Time	Set date, time and display format. Will be used for the Browse function.
Color Scheme	Change color theme on display (bright / dark).
Information	View current firmware version and serial number for instrument.
Save	Turn saving of results on or off in Viscosity Measurement .
Lock Settings	Lock current settings against unintentional changes.
Table Settings	Select table format for measuring screen and the number of readings per second / minute.
Change Cone	View list of existing cones. Select the cone to be used for measuring.
Averaging	Off: The system records a measured value every 0,5 seconds. The result may look unstable. On: The last "n" measured values are averaged. The result may look more stable.
Torque Range	Low: Max. torque value = 0.797 mNm. High: Max. torque value = 18.1 mNm.

NOTICE

Configuration for **Units** will be used as default and can also be changed directly during measurements.


6 Viscosity Measurement

The measurement takes place in two steps - heating up (or cooling down) and measuring:

1. Tap the green cone icon lower the viscometer head the 1st time.
2. Remove any sample completely from the plate; confirm the message.
3. Viscometer head moves down, the system regulates the temperature.
4. When target temperature is reached, confirm the message to raise the cone.
5. Apply the sample to the plate.
6. Tap the blue cone icon to lower the viscometer head the 2nd time.
7. Play button becomes green; tap to start the measurement.


Heating up to high temperatures may take some time.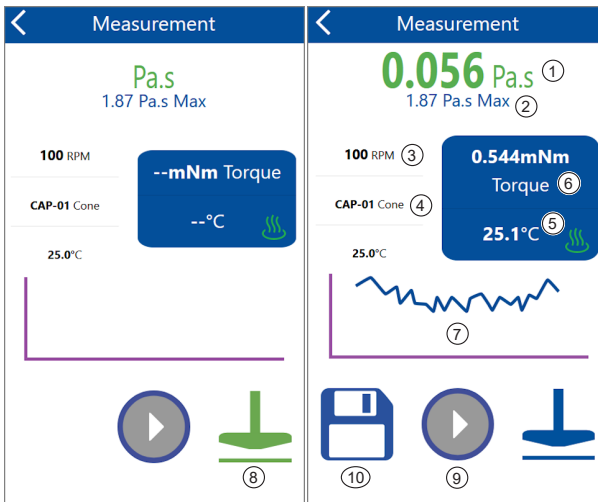

1	Viscosity	Value is taken every 0.5 seconds; tap to change unit.
2	Maximum	The max. value depends on cone, speed and torque setting.
3	Speed	Current cone speed in rotations per minute, tap to change.
4	Cone	Identifier for current cone, tap to change.
5	Temp.	Current temperature of sample: Green: OK. Red: Heating up. Tap to change unit.
6	Torque	Current value mNm or in percent of max. value; tap to change; change speed or cone or torque setting, if < 15% or > 95% (invalid measurement).
7	Graph	Shows a live updating graph of viscosity over time.
8	Lower	Green: Tap to regulate temperature. Blue: Tap to measure.
9	Start	Green: Start measurement. Gray: Inactive.
10	Disk	Save results in system memory; appears after stopping measurement.

Touch areas of the screen with measured data to access corresponding setting.

Swipe left and right to access the other screens.

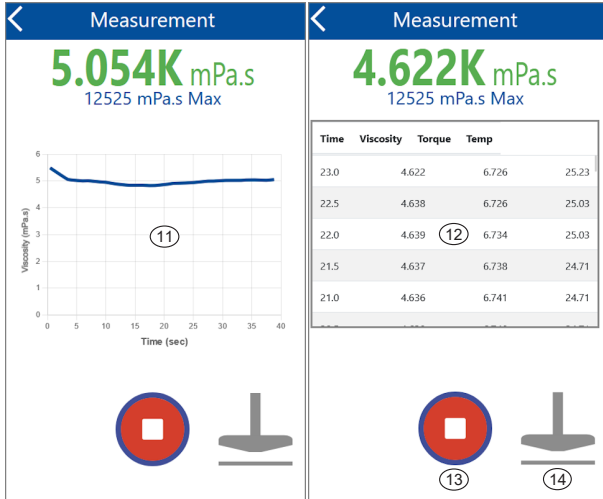
- 11 Graph Shows measurement data in graphical form.
- 12 Table Shows measurement data in table form.
- 13 Stop Red: End measurement and save results.
- 14 Lower Gray: Inactive.

7 Browse Measurements

Tap the browse icon in the main menu to manage saved measurement data. The final value for a measurement is displayed here. Export to USB to see all viscosity values measured over time.

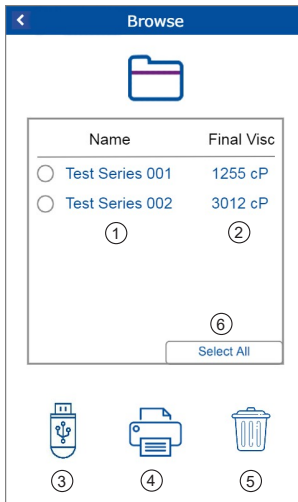1	File Name	Identify a test series by its name, tap to rename.
2	Final Visc	Gives the final viscosity value for this test series.
3	Write USB	Write selected test series to USB flash drive.
4	Print USB	Print selected test series to USB printer.
5	Delete	Delete selected test series from memory.
6	Select All	Select / deselect all rows / test series at once.

Active options are dark blue. Inactive options are light blue.

8 Technical Data

Instrument

Catalog number	5210: byko-visc CP L (Low temperature) 5211: byko-visc CP H (High temperature)
Temperature range	byko-visc CP L: 5° C to 75° C byko-visc CP H: 50° C to 235° C
Torque range	Low: Maximum = 0.797 mNm High: Maximum = 18.1 mNm
Shear range	10 to 13,000 sec ⁻¹
Speed range	5 to 1,000 RPM
No. of speeds	Freely selectable within speed range
Repeatability	0.2%
Accuracy	2% of full-scale range
Preset memory	6 memory locations for measurement settings presets
Measurement features	Time to Stop, Multi-step Ramp, QC Limits, Curve Fitting, Multi-point Tests
Display	Capacitive color display, 480 x 800 pixels
Languages	English, German, French, Italian, Spanish, Russian, Japanese, Chinese

General

Temperature range	10° C to 40° C (50° F to 104° F) for operation 0° C to 60° C (32° F to 140° F) for storage
Relative humidity	Up to 85% non-condensing at 35° C (95° F)
Operation altitude	Up to 2000 m (6561 ft)
Dimensions (LxWxH)	300 x 170 x 430 mm (11.5 x 6.5 x 17 in)
Weight	6.5 kg (14.3 lbs)
Interface	2x USB Type A 1x USB Type C 1x Ethernet
Power supply instrument	byko-visc CP L models: 24 V  , max. 3.75 A byko-visc CP H models: 24 V  , max. 5 A
External power supply	byko-visc CP L models: Input: 100 - 240 V  ; 50 - 60 Hz ; max. 1.3 A Output: 24 V  ; max. 3.75 A ; max. 90 W byko-visc CP H models: Input: 100 - 240 V  ; 50 - 60 Hz ; max. 1.4 A Output: 24 V  ; max. 5 A ; max. 120 W

Technical data are subject to change without notice.

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BYK-Gardner GmbH