## byko-swing

Short Instructions

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Instructions rapides

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Instrucciones resumidas

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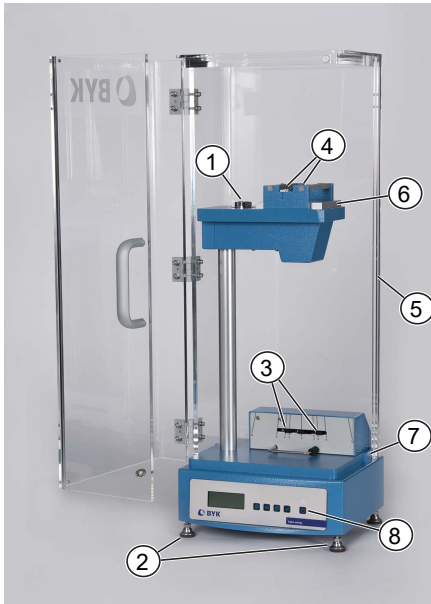
# Table of Contents

|                |    |
|----------------|----|
| English .....  | 4  |
| Deutsch.....   | 6  |
| Français.....  | 8  |
| Italiano ..... | 10 |
| Español .....  | 12 |
| 中文 .....       | 14 |
| 日本語 .....      | 16 |

# 1 System Description

The system is supplied fully assembled - including protective cover, pendulum(s), glass calibration standard and power supply.

|                                                            |                               |
|------------------------------------------------------------|-------------------------------|
| 1 Bubble level                                             | 2 Leveling feet               |
| 3 Light barriers                                           | 4 Pins for pendulum           |
| 5 Protective cover                                         | 6 Automatic lifting table     |
| 7 Door sensor to ensure cover is closed during measurement | 8 Power switch and status LED |

1. Place the instrument on a vibration free surface.
2. Connect the external power supply, socket on back site.
3. Balance the bubble level (1) properly with the levelling feet (2).
4. Place the pendulum into the tapered pins (4) of the lifting table (6), **arrow pointing to the front.**
5. Close the door of the protective cover, double check bubble level.


## NOTICE

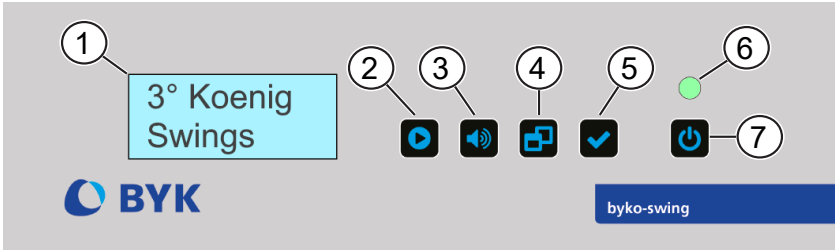The device is heavy - transports are to be done with the help of a second person. Retain the packaging in case the instrument needs to be shipped.

## 2 Main Menu and Operation

To ensure accurate measurements, clean the housing with isopropanol before use. For more information, please refer to the Operating Manual.

### Operating the Instrument

- Power on the instrument (7).
- Place sample on automatic lifting table.
- Clean the balls of the pendulum with isopropanol.
- Select appropriate method with button **Modus** (4).
- Confirm with **Enter** (5), status LED (6) lights up green for *Ready*.
- Optional: Adjust duration of acoustic end signal with button **Sound** (3) and confirm with **Enter** (5).
- Press button **Start/Stop** (2), the LED lights up red for *Busy*.

The following four measurement modus are available by pressing the button **Modus** (4).

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The swing time for "König" is 1.4 seconds, for "Persoz" it is 1.0 second.

### Testing with Glass Calibration Standard

It is recommended to check the accuracy with the included glass calibration standard **once a week**:

1. Clean the glass calibration standard by soft, lint-free cloth with isopropanol. Use always a new cloth for cleaning.
2. Place glass standard on automatic lifting table and start measurement.
3. If the values cannot be obtained, reclean the glass panel and the balls of the pendulum, recheck that the glass panel is levelled, and retest.

|                            | Swings (n) | Time (s)   |
|----------------------------|------------|------------|
| Pendulum hardness - König  | 172 - 185  | 250 +/- 10 |
| Pendulum hardness - Persoz | 415 - 445  | 430 +/- 10 |

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**BYK-Gardner GmbH**